ZeGlobalTox, a new way for global toxicity assessment

Carles Cornet, Rafael Miñana, Simone Calzolari, Davide D'Amico and Javier Terriente. ZeClinics, PRBB, Barcelona, Spain, www.zeclinics.com.










Zebrafish exploitation for high-throughput drug screening is becoming important for the pharmaceutical industry to assess toxicity and efficacy of novel drugs. Zebrafish has, from early developmental stages, fully functional organs from a physiological point of view. Thus, drug-induced toxicity is easily detected and evaluated, bridging the gap between preclinical in vitro and in vivo models in a fast and cost-effective manner.

ZeClinics has developed ZeGlobalTox, an innovative tool that integrates in vivo toxicity assessment for the brain, heart and liver in the same animal, strongly impacting the 3Rs principles. This assay *Reduces*, by up to a third, the number of animals required to assess toxicity in these organ. The new physiological parameters included in the ZeGlobalTox Refines the drug toxicity assay. The high predictivity of the ZeGlobalTox (Specificity: 85% and Sensitivity: 76%) allows to *Replace* the use of other classical species, such as rodents or other mammals.

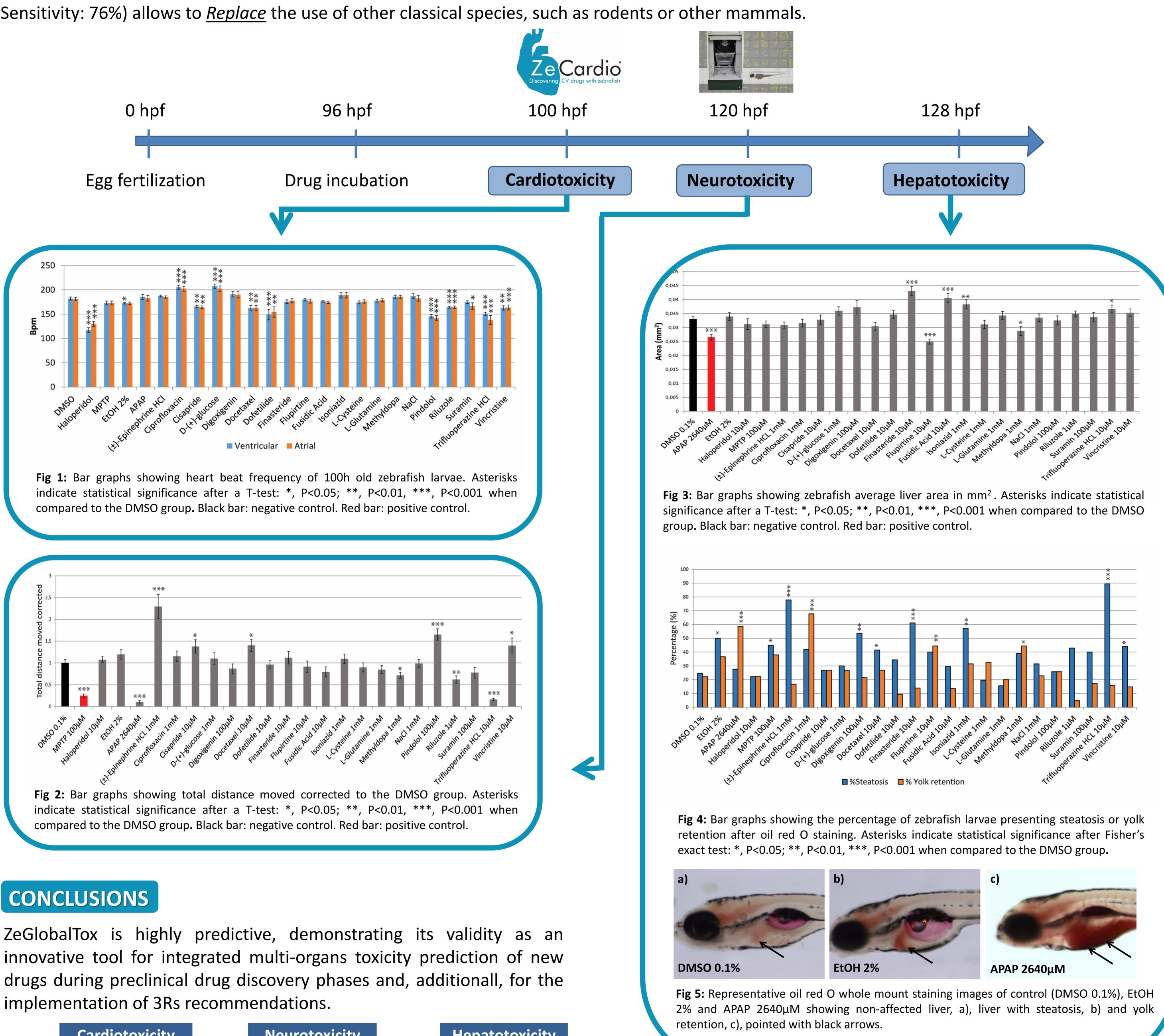

implementation of 3Rs recommendations.

	Cardio		oxicity			Neurot	oxicity			Hepatotoxicity	
		Human				Hum	nan			Human	
		-	+			-	+			-	+
Zf	-	TN: 12	FN: 2		- TN: 13	FN: 2		-	TN: 17	FN: 2	
	+	FP: 3	TP: 8	Zf	+	FP: 2	TP: 8	Zf	f +	FP: 2	TP: 4

Zf: zebrafish, **TN:** true negative, **TP:** true positive, **FN:** false negative, **FP:** false positive.

	Specificity TN/(TN+FP)	Sensitivity TP/(TP+FN)	PPV TP/(TP+FP)	NPV TN/(TN+FN)
Cardiotoxicity	80%	80%	73%	86%
Neurotoxicity	87%	80%	80%	87%
Hepatotoxicity	89%	67%	67%	89%
ZeGlobalTox	85%	76%	73%	87%

Zf: zebrafish, **PPV**: positive predictive value, **NPV**: negative predictive value, **TN:** true negative, **TP**: true positive, **FN**: false negative, **FP**: false positive.

ACKNOWLEDGEMENT & FUNDING

The ZeGlobalTox project was run in collaboration with the Pivot Park Screening Center, Oss, The Netherlands, which provided the set of tested compounds and Noldus Information Technology, Wageningen, The Netherlands, contributing by providing the DanioVision TM for the behavioral analysis.

Financial supports have been received by the National Centre for the Replacement, Refinement and Reduction of Animals in Research (NC3Rs) as an innovative CRACK IT Solution – Agreement of June 1st 2015 - and by the Spanish Ministry of Economy and Competitiveness though a Torres Quevedo Fellowship (PTQ-14-07351) to SC and an Industrial Doctorate Fellowship (DI-14-06969) to CC.





